


Anesthesia Mask Connections

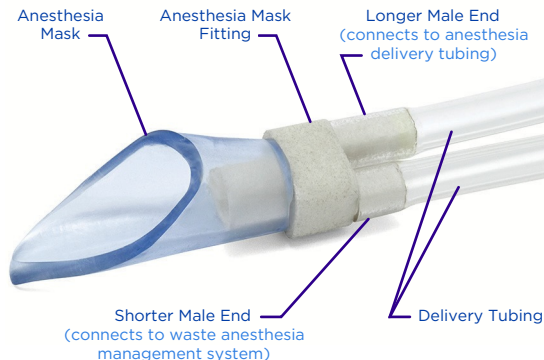

Anesthesia Mask

Each kit includes one anesthesia mask already attached to the fitting and tubing.

NOTE: The longer male end on the back of the mask fitting identifies the anesthetic delivery tubing which should always be connected to the vaporizer. The shorter male end should be connected to your waste anesthesia gas management system.

Tips for a Good Fit

- The mask can be used in any orientation; simply rotate the mask on the white anesthesia mask fitting.
- Ensure the animal's nose and mouth are covered.
- The mask should fit snugly around the animal's muzzle without making contact with the eyes.

Vaporizer Connections

Traditional VetFlo™

Includes:

36" delivery tubing

1 set (2 pieces) 1/4" tube adapters

1 set (2 pieces) 3/8" tube adapters

Select the appropriate sized adapter to fit to your vaporizer.

SomnoSuite®

Includes:

18" delivery tubing

Male Luer connectors **(2)**